
WCAC AGILENT 1100 HPLC USAGE RULES

10AUG2014

THE AGILENT HPLC RATE: \$1:00/30MIN RESERVATION TIME

SIGN IN: A) NAME B) NETID, C) LAB PI, D) CONTACT INFO IF YOU ARE NOT PRESENT DURING THE ENTIRE RUN E) NUMBER OF RUNS AND TOTAL ON TIME.

AGILENT 1100 HPLC USE:

1. Reserve at <http://scheduler.yale.edu>, for cancellations email WCAC staff.
2. Columns and solvents are the user's responsibility.
3. **Use the instrument only after training by WCAC Staff.**
4. Prime pumps/purge with your own solvent, if necessary.
5. At the end of your run, purge the pumps with 50% Isopropanol/water to preserve the seals.
6. Properly dispose of the flow through/waste effluent in YOUR parent lab.
7. Shutdown the system when done.

Please alert staff of any instrument faults/operating trouble (e.g., leaks, flow, pressure).